

## Overview

It is often useful to have a command-line utility that simply triggers the TG6 at regular intervals. The TG6\_SW2.EXE program was written to trigger up to six TG6 channels and associated SW2s. You can use also use it without any SW2s.

## Usage

To use the program, first type the program name at the DOS prompt without any arguments. The program prints out a list of the arguments it requires.

The program automatically rounds values to the resolution of timing that has been selected and range-checks for some (but not all) errors that can occur.

Running TG6\_SW2 with one argument turns off the TG6 (e.g. "TG6\_SW2 off").