

Medusa4Z BioAmp

Hardware Reference



Updated 2025-02-27

© 2016-2025 Tucker-Davis Technologies, Inc. (TDT). All rights reserved.

Tucker-Davis Technologies
11930 Research Circle
Alachua, FL 32615 USA
Phone: +1.386.462.9622
Fax: +1.386.462.5365

Notices

The information contained in this document is provided "as is," and is subject to being changed, without notice. TDT shall not be liable for errors or damages in connection with the furnishing, use, or performance of this document or of any information contained herein.

The latest versions of TDT documents are always online at <https://www.tdt.com/docs/>

Table of Contents

Medusa4Z BioAmp

Medusa4Z Overview	4
Medusa4Z Features	5
Analog Input Channels	5
Electrode Connectors	6
Impedance Checking	6
LCD Screen	6
Base Station Connector - Optical	7
Power	8
Power Requirements	8
Medusa4Z Technical Specifications	9
Input Referred Noise	10
Sample Delay	10

Medusa4Z BioAmp



Medusa4Z Overview

The Medusa4Z is a four-channel, referential, low-noise digital bioamp targeted for use in low-channel-count, low-impedance neurophysiology recordings, such as ABR, EEG, EMG, or EKG signals. The bioamp digitizes one or four channels at acquisition rates of approximately 3 kHz, 6 kHz, 12 kHz, or 25 kHz, and the amplified digital signal is sent to the base station via a noiseless fiber optic connector.

Connections to the subject (four channels, reference, and ground) are made with six standard 1.5 mm touchproof safety connectors.

- It is powered by a 5 AmpHr Li-Poly battery that provides up to 30 hours of continuous data acquisition in single-channel mode and up to 24 hours in four-channel mode.
- It includes impedance testing functionality for each of the four recording channels as well as the reference electrode.
- The user interface has 8 buttons and a 16 x 2 character LCD display screen. It lets you select performance characteristics of the preamplifier and shows battery and signal level.
- A complete charge from a dead battery can be achieved in under 8 hours with the included charger.

Medusa4Z Features

The Medusa4Z has the option to be run in single-channel mode or four-channel mode. When in single channel mode, channels 2-4 are powered down and isolated from the reference line. Running in single-channel mode will provide slightly better performance and longer battery life.

There are selectable Sampling Frequency options of ~3 kHz, ~6 kHz, ~12 kHz, or system rate (maximum ~25 kHz). Users will find that using the lowest practical rate improves noise performance.

When testing the impedance of the probes, users can select frequencies of 100 Hz, 300 Hz, 500 Hz, or 1 kHz.

A user-selectable digital highpass filter has adjustable options that are: Off, 1 Hz, and 3 Hz. The hardware has a fixed analog highpass at 0.3 Hz.

An Auto Shutdown feature has three modes: Never, 1 hour, 3 hours, or 5 hours. If no buttons are pressed after this interval, then the device will automatically power down. If no optics are detected, the Medusa4Z will also shutdown. Before shutting down, the blue POWER LED will blink fast to indicate that time has expired. Pressing any key will reset the counter.

Analog Input Channels

The Medusa4Z Preamplifier acquires signals with 16-bit PCM ADCs. The signals are oversampled before conversion to remove aliasing of high frequency RF signals. This allows users to acquire at lower sampling rates (down to 3 kHz) with improved signal-to-noise quality in that range and no aliasing. The Medusa4Z can also acquire data in High-Resolution mode which uses 24-bit PCM ADCs. This is controlled by the base station (RZ, RX) to which the Medusa4Z is connected. The change occurs by modifying the underlying circuit running on the base station in Synapse or BioSigRZ. This is handled by default in [Synapse v92](#) and above, and [BioSigRZ v5.7.5](#) and above. In TDT Drivers v92 and above there is a RPvdsEx circuit macro available in: `C:\TDT\RPvdsEx\Macros\Device\Medusa4Z_Bioamp`.

Electrode Connectors

Plug the electrodes directly into the standard 1.5 mm touchproof connectors on the side of the Medusa4Z.



Impedance Checking

The impedance checker on the Medusa4Z provides a simple check of the channel impedance relative to ground. To check the impedance level, press the button next to the channel indicator (e.g. CHAN-1). The impedance between channel and ground will be shown on the screen. The REF impedance button checks the impedance between the reference and the ground. Impedance checking frequency options are 100 Hz, 300 Hz, 500 Hz, and 1 kHz. These are settable on the user interface.

LCD Screen



The LCD screen shows the current sampling rate, battery level, and signal level indication on the active channels.

Use the SETUP button to cycle through available settings and use the OPTION button to change the currently selected setting. Settings include:

Active Channels: Single or Four Channel

Sample Frequency: ~3 kHz, ~6 kHz, ~12 kHz, System Rate (up to ~25 kHz). Match this to the desired frequency range of your acquired signal (maximum frequency captured will be ~45% of selected sample frequency). Note that lower sampling rates will have improved signal-to-noise, but they will also have longer sample delays. Please refer to the See Medusa4Z PreAmp Technical Specifications. for more information.

Probe Frequency: 100 Hz, 300 Hz, 500 Hz, 1 kHz. Choose the default frequency used for the impedance checker.

Highpass Filt: Off, 1 Hz, 3 Hz. Set the highpass filter cutoff frequency on the input signals. 'Off' is a 0.1 Hz filter. Note that there is a fixed 0.3 Hz highpass analog filter on the inputs.


Auto Shutdown: Never, 1 hr, 3 hr, 5 hr. After a period of inactivity the Medusa4Z will shut itself down to preserve power. This prevents dead batteries from accidentally leaving the unit on overnight.

Base Station Connector - Optical



Use the provided fiber optic pair (black cable, white connectors) to connect the OPTICAL port on the preamplifier to the optical input port on the base station.

The duplex fiber optic cable has identical one-piece connectors at each end. There is a V-shaped groove on one side of the connector and a raised rectangle on the other. As shown in the image below, plug the connector into the RZ6 port with the raised rectangle side up. Plug the connector into the preamplifier with the V-shaped groove up.

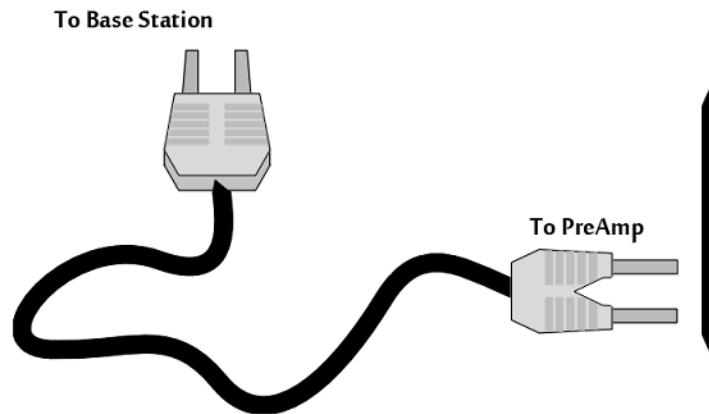

Medusa4Z Connection Diagram


Power

The POWER button on the face of the Medusa4Z powers it on. Press and hold the button to power it off. Press the power button momentarily to go to the home screen from anywhere or to reset the power down timer.

Do not use the recessed slide power switch (ON OFF (RESET)) on the side of the device unless you need to reset all settings to factory default, or if you are storing the Medusa4Z for an extended period of time and want to completely disconnect the battery. The OPTICAL fiber connector at the right will be illuminated when the amplifier is on.

Power Requirements

The lithium battery charges in four hours. Keeping the battery charger connected to the amplifier does not affect the battery life. However, the charger will significantly increase the noise of the system if it is plugged in while an experiment is running. A 12 Volt battery charger is included with the amplifier. The supplied charger tip is center negative. However, tip polarity can be positive or negative if the charger needs to be replaced.

 **Note**

If the Medusa4Z is powered on while charging, the charge circuit will halt when the battery is full. It will resume charging when the battery runs down to ~90%.

The lithium battery supplied with the system cannot be removed. If battery life longer than 30 hours is required, an external battery pack can be connected to the voltage inputs of the charger. TDT recommends a 6 V (minimum) to 12 Volt (maximum) battery, such as lead acid batteries used for motorized wheel chairs. Contact TDT for more information.

Medusa4Z Technical Specifications

A/D	Up to 4 channels Standard Mode: 16-bit PCM High-Resolution Mode: 24-bit PCM
Sample Rate	~3 kHz, ~6 kHz, ~12 kHz, or ~25 kHz
Maximum Voltage In	±10 mV
Gain	100x
Frequency Response	3 dB 0.3 Hz - 0.45*Fs (with no additional high-pass filter selected)
Highpass Filter	0.3 Hz, 1 Hz, or 3 Hz
Anti-Aliasing Filtering	45% sampling rate
S/N (typical)	80 dB
Input Impedance	100 kOhm
Battery	Li-Poly Battery 5000 mAh single-channel mode: 30 hour runtime four-channel mode: 24 hour runtime 1000 cycles of charging, not removable by user
Charger	6-16 V DC, >1 A, center-positive or center-negative
Fiber Optic Cable	5 meters standard, maximum cable length 12 meters

Input Referred Noise

Depends on Medusa4Z sample rate and electrode impedance. All units in μVrms .

Medusa4Z rate	shorted	3 kOhm	10 kOhm
3 kHz	0.15	0.3	0.5
6 kHz	0.25	0.4	0.7
12 kHz	0.3	0.6	1.0
25 kHz	0.4	0.9	1.4

Sample Delay

Depends on Medusa4Z and RZ processor sample rates. All units in samples.

Medusa4Z rate	RZ @ 25 kHz	RZ @ 12 kHz	RZ @ 6 kHz
25 kHz	24	N/A	N/A
12 kHz	42	24	N/A
6 kHz	76	42	24
3 kHz	146	76	42