

iL24 Digital Logic Interface

Hardware Reference



Updated 2024-02-05

© 2016-2024 Tucker-Davis Technologies, Inc. (TDT). All rights reserved.

Tucker-Davis Technologies
11930 Research Circle
Alachua, FL 32615 USA
Phone: +1.386.462.9622
Fax: +1.386.462.5365

Notices

The information contained in this document is provided "as is," and is subject to being changed, without notice. TDT shall not be liable for errors or damages in connection with the furnishing, use, or performance of this document or of any information contained herein.

The latest versions of TDT documents are always online at <https://www.tdt.com/docs/>

Table of Contents

iL24 Digital Logic Interface

DB25 Pinout

iL24 Digital Logic Interface




The iL24 module can communicate with external behavioral components using 24 bits of 5 V or 3.3 V TTL logic signals. There are 12 input bits and 12 output bits. Each direction has 4 addressable bits and 1 addressable byte.

- There are four bit-wise inputs and four bit-wise outputs available with BNC connectors and status light for each
- The DB25 connector has access to all 24 addressable bits
- Two rows of 8 status LEDs on the front panel show the state of the word input and output bits
- The front panel switch toggles between +3.3 V and +5 V logic for all 24 bits of I/O on the iL24
- Each bit can source up to 6 mA maximum current

For information on software control of the iL24, see the [Synapse Manual](#).

DB25 Pinout

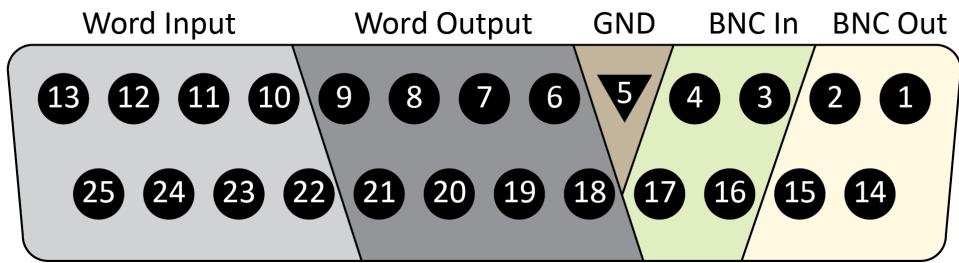

iL24 pinout, looking into the connector

Pin	Name	Description	Pin	Name	Description
1	B01	BNC Output	14	B02	BNC Output
2	B03		15	B04	
3	B11	BNC Input	16	B12	BNC Input
4	B13		17	B14	
5	GND	Ground	18	D00	Word Output
6	D01	Word Output	19	D02	
7	D03		20	D04	
8	D05		21	D06	
9	D07		22	D10	Word Input
10	D11	Word Input	23	D12	
11	D13		24	D14	
12	D15		25	D16	
13	D17				