

# iL24 Digital Logic Interface

---

Hardware Reference



Updated 2025-02-27

© 2016-2025 Tucker-Davis Technologies, Inc. (TDT). All rights reserved.

Tucker-Davis Technologies  
11930 Research Circle  
Alachua, FL 32615 USA  
Phone: +1.386.462.9622  
Fax: +1.386.462.5365

### **Notices**

The information contained in this document is provided "as is," and is subject to being changed, without notice. TDT shall not be liable for errors or damages in connection with the furnishing, use, or performance of this document or of any information contained herein.

The latest versions of TDT documents are always online at <https://www.tdt.com/docs/>

# Table of Contents

---

## **iL24 Digital Logic Interface**

iL24m Manifold	4
DB25 Pinout	5

# iL24 Digital Logic Interface

---



The iL24 module can communicate with external behavioral components using 24 bits of 5 V or 3.3 V TTL logic signals. There are 12 input bits and 12 output bits. Each direction has 4 addressable bits and 1 addressable byte.


- There are four bit-wise inputs and four bit-wise outputs available with BNC connectors and status light for each
- The DB25 connector has access to all 24 addressable bits
- Two rows of 8 status LEDs on the front panel show the state of the word input and output bits
- The front panel switch toggles between +3.3 V and +5 V logic for all 24 bits of I/O on the iL24
- Each bit can source up to 6 mA maximum current

For information on software control of the iL24, see the [Synapse Manual](#).


## iL24m Manifold

---

The iL24 comes with a manifold of BNC connectors for easier access to the Word Input and Word Output bits.

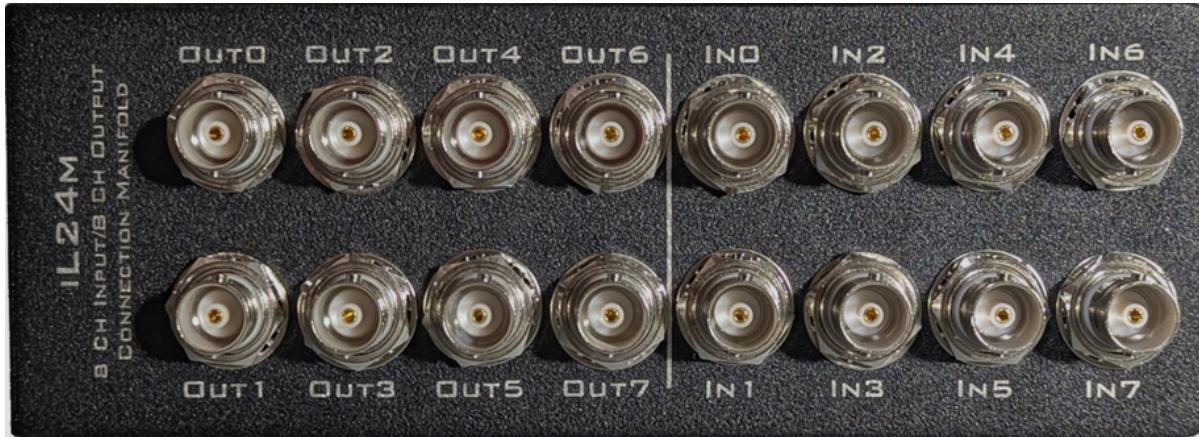

## DB25 Pinout

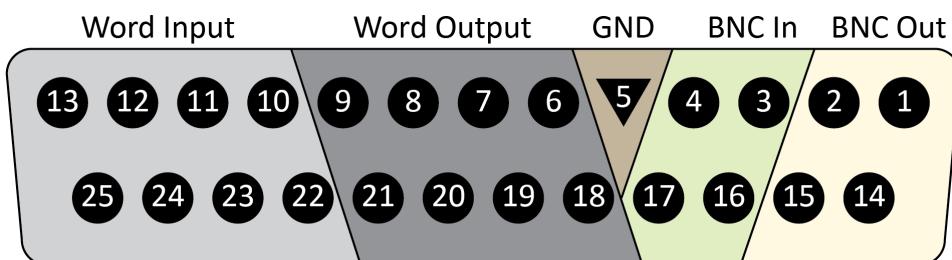

iL24 pinout, looking into the connector

Pin	Name	Description	Pin	Name	Description
1	BO1	BNC Output	14	BO2	BNC Output
2	BO3		15	BO4	
3	BI1	BNC Input	16	BI2	BNC Input
4	BI3		17	BI4	
5	GND	Ground	18	DO0	Word Output
6	DO1	Word Output	19	DO2	
7	DO3		20	DO4	
8	DO5		21	DO6	
9	DO7		22	DI0	Word Input
10	DI1	Word Input	23	DI2	
11	DI3		24	DI4	
12	DI5		25	DI6	
13	DI7				